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General Regulations of the laboratories of experimental psychology (Experimental Psychology Lab) , CIMeC - Fedrigotti Palace, Rovereto

December 2013

Additional Rules for the use of laboratory and instrumentation for the study of EYE MOVEMENTS

The supplement of regulation provided for the use of the equipment for the study of eye movements is due to the fact that the devices are sophisticated and expensive. The equipment must be handled with particular care. For its use some competence is needed.

The eye-tracking system (EyeLink 1000 SR -research) requires a system of monitoring / recording "host PC" and is used in connection to a PC for the generation of stimuli (computer display) . The communication between the two computers is via Ethernet and you need to change the network settings of the PC stimulation and install a special software package.

The various systems are equipped with an identification code associated with a file that must be added to the management software of the host, so for the system to work with a given host the file associated with the device that you want use must be installed. Although very similar, the EyeLink systems in use have small differences that make it necessary to use a management software for recording, or specific (host) computers. This does not influence the programming of the experiments, which remains the same for the various systems.

In some laboratories there are specific workstations ("Display PC" connected to a "host PC" , to which you only have to connect the optics) to be used for experiments in which you want to monitor eye movements. The various optical and portable systems are stored in a common space; information about available resources, their location and how to use the various systems are provided in the common space.

For the Tobii (Tobii T120), there is a single computer for recording of eye movements and the administration of the stimuli; the software, Tobii Studio, is installed on a laptop, in order to make the system more easily portable. There are no default locations in which to use the Tobii.

- To start a new study that requires the use of one or more systems for the control of eye movements or if there is the need to modify the operation, the arrangement or positioning of one of the Eyetracking systems, you must request a preliminary meeting with the technician in charge of the laboratories.

CiMeC

Palazzo Fedrigotti
corso Bettini, 31 - 38068 Rovereto, Trento (Italy)
ph. +39 0464 808615, fax +39 0464 808690
e-mail: cimec@unitn.it
<http://www.cimec.unitn.it>

- Handle with care optics and connection cables: even a small damage can cause time and high intervention costs. The optics and various parts of the device should not be removed unless previously agreed with the technician.
- Do not modify the software or initialization files (INI files) on the recording PC (host) except in accordance with presence of the technician. Any accidental modification of the software or recording settings must be communicated promptly to the technical and other users.
- To reduce the wear of the devices (especially the connectors), and the possibility of accidental damage, it is appropriate to limit frequent moves and connections/disconnections. For this reason it is advisable to condensate as much as possible the data collection sessions and to agree on the use of equipment with other users.
- If the device is not used for a long period, it must be stored in the appropriate space. When the device is used regularly it is preferable to leave it connected to the appropriate workstation to avoid repeated connections / disconnections.

Safety Rules

The infrared radiation emitted by infrared illuminators installed on the devices can be harmful to the eyes. For this reason, it is important to avoid staring at the LEDs from a close distance.